## Compact High Performance

# Sirius L Single Tube Luminometer

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- Reporter gene assays, including dual-reporter assays
- Rapid detection of mycoplasma contamination in cell cultures: Lonza MycoAlert<sup>®</sup>
- ATP assays
- Cell proliferation, cytotoxicity and biomass assays
- Apoptosis assays
- Cellular luminescence/reactive oxygen species
- Environmental toxicity and mutagenicity assays

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- 1.5 ml, 2 ml microfuge tubes
- 75×12 mm, 55×12 mm tubes
- 35 mm petridish,

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# Compact High Performance Sirius L Single Tube Luminometer

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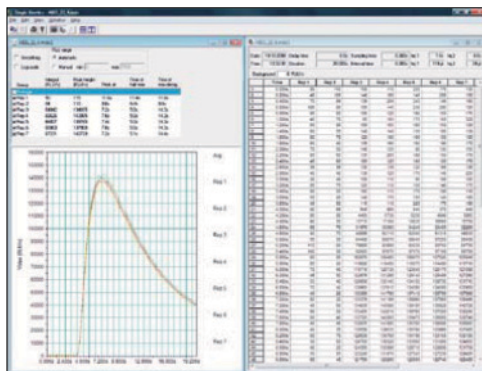
그림 5. Available protocols for every application, Kinetic evaluation, data table (right), curve parameters and graphic presentation (left)

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### Luminometer

Detector	Photomultiplier tube with bialkali cathode, effective spectral range 300–630 nm, operated in photon counting mode.
Sensitivity	Less than 0.5 zmol firefly luciferase, 1 amol ATP in a HS ATP assay.
Dynamic Range	6 decades.
Sample Format	Tubes up to 12 mm diameter and 75 mm length, microfuge tubes 1.5 ml and 2 ml, 35 mm culture dishes, liquid scintillation vials up to 20 ml. Injection operation with tubes 12 mmx75 or 55 mm and microfuge tubes of 1.5 and 2 ml volume.
Measurement Chamber	Retractable drawer with interchangeable reflectors and sample adapter.
Injector	0, 1, 2 (upgrade possible).
Injector Volume	20-500 $\mu$ l adjustable in steps of 1 $\mu$ l.
Tubing	Chemically inert PTFE tubing and connections (PTFE or KEL-F).
Interfaces	RS 232 or USB interface for PC connection.
Dimensions	W 349 mm, D 150 mm, H 241 mm (without injection tubings).
Weight	approx. 3,6 kg (5,4 kg with 2 injectors).
Power Requirement	Desktop power supply and power cord. 100-240 V AC, 24 V/2,5 A DC.
Environment	Storage 0-40° C up to 80% humidity (30° C), non condensing. Operation 10-35° C up to 70% humidity.

### PC Software

PC Software	Microsoft Windows® application for control, measurement and evaluation.
Required Hardware	Microsoft Windows® compatible PC, Pentium like processor, RS 232 or USB port.
Operating System	Windows® 2000, XP, Vista, Windows® 7.
Additional Software	Microsoft Excel™ 2000 or higher (optional).
Standard Configuration	Protocol Manager, Quick Measurement.
Additionally Available	Single Assay, Dual Assay, Single Kinetics, Multiple Kinetics, Cut-off Assay.