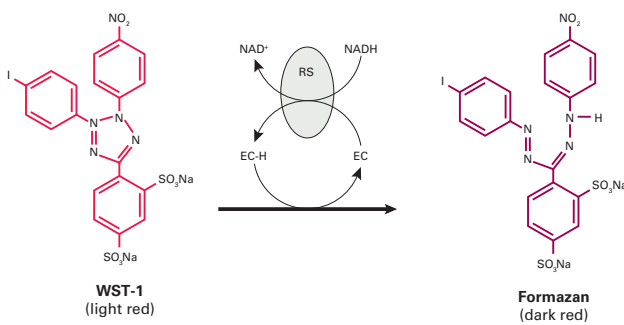
# Utilizing the Premixed WST-1 Cell Proliferation Reagent to Avoid Off-Target Effects of RNAi

Peter W. Roelvink\* and David A. Suh\*  
 \* Currently with Clontech Laboratories, Inc.

Work performed while at Benitec Ltd., Hawthorne East, Australia

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**그림 1. Chemical structure and cleavage of the tetrazolium salt WST-1 to formazan.** Once in the interior of active cells, WST-1 is reduced by mitochondrial succinate-tetrazolium reductase to produce the colorimetrically differentiated formazan product. EC = electron coupling reagent. RS = mitochondrial succinate-tetrazolium reductase system

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
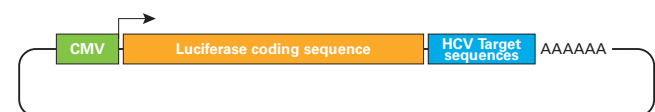
**그림 2. Luciferase reporter system for assaying RNAi activity.** The reporter was constructed by fusing HCV sequences (~100 nucleotides of contiguous viral sequence) downstream of the luciferase reading frame but upstream of the poly A sequence.

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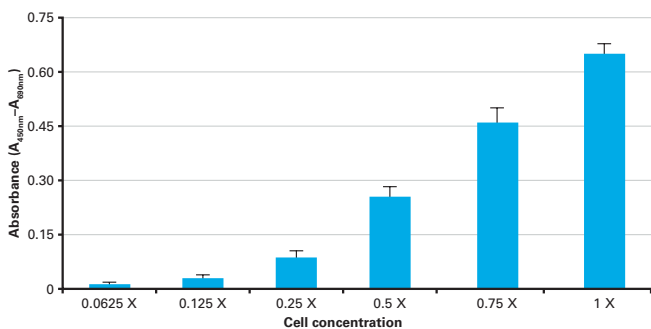
**그림 3. Ability of WST-1 reagent to detect differences in proliferating cells.** Huh-7 cells were plated at differential densities and allowed to grow for an additional 72 hr prior to performing the WST-1 assay. Measurements are the averages of n = 4 independent samples ± standard deviation.

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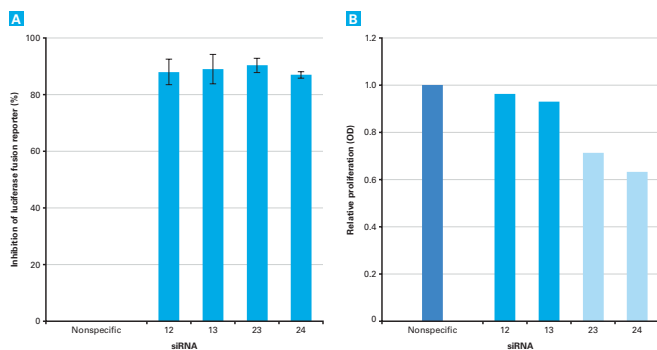**그림 4. Assessment of nonspecific siRNA toxicity using WST-1.** Panel A. Inhibition of specific luciferase-HCV reporter plasmid activity in Huh-7 cells 72 hr after cotransfection with the cognate siRNA (n = 4 independent transfections). The data were normalized for transfection efficiency by inclusion of a second expression plasmid encoding Renilla luciferase in the transfection mixture. Inhibitory activity was calculated based on comparisons of parallel transfections that utilized a nonspecific control siRNA. The data are represented as the mean inhibition of luciferase activity relative to the nonspecific siRNA control ± standard deviation. Panel B. Assessment of the effects of the siRNA on cellular proliferation by the WST-1 assay. Absorbance readings were normalized to cells that had been transfected with the nonspecific control siRNA (n = 4 independent transfections).

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Code	제품명	용량
630118	Premixed WST-1 Cell Proliferation Reagent	2,500 rxns
631726	Ready-To-Glow Secreted Luciferase Reporter Assay	100 회
631727	Ready-To-Glow Secreted Luciferase Reporter Assay	500 회
631728	Ready-To-Glow Secreted Luciferase Reporter Assay	1,000 회
631729	Ready-To-Glow Secreted Luciferase pMetLuc Vector Kit	1 kit

License Notice [K24]

**References :**

- Birmingham, A. *et al.* (2006) *Nat. Methods* 3(3):199-204.
- Semizarov, D. *et al.* (2003) *Proc. Natl. Acad. Sci. USA* 100(11):6347-6352.
- Jackson, A.L. *et al.* (2003) *Nat. Biotechnol.* 21(6):635-637.
- Slater, T.F. *et al.* (1963) *Biochim. Biophys. Acta* 77:383-393.
- Kittler, R. *et al.* (2004) *Nature* 432(720):1036-1040.